

Reference Review  
Spore News

## Biological Indicators for Vaporized Hydrogen Peroxide Isolator Decontamination: Characteristics, Uses and Challenges

The process of validating an isolator decontamination process was discussed in a recent Mesa white paper ([Vaporized Hydrogen Peroxide Isolator Decontamination in a World of Uncertainty](#)), and therein, the selection and qualification of biological indicators (BIs) was covered in moderate detail. This Spore News Reference Review document is entirely dedicated to vapor hydrogen peroxide ( $\text{VH}_2\text{O}_2$ ) BIs including necessary physical characteristics of the BI unit, BI types, proper BI placement, and challenges encountered.

It is recognized that  $\text{VH}_2\text{O}_2$  has poor penetration abilities; however, well-designed processes have demonstrated that it can be an effective agent for surface decontamination, provided the methods/equipment control the process within the parameters empirically proven to be effectively lethal.

One key prerequisite to the decontamination process is a cleaning step. The purpose of cleaning the isolator is to remove materials that may adversely affect the product to be produced in the system. These materials can be general debris along with microbial contamination. Even the most robust cleaning process cannot remove all contaminating bioburden; however, microorganisms that are still present after the cleaning are made more vulnerable to the decontamination process. This is accomplished by removing any large aggregates of microorganism and the materials that may shield the remaining organisms during the decontamination process. Likewise, BIs used to monitor  $\text{VH}_2\text{O}_2$  processes must present the spores in a similar fashion as the organisms remaining in the isolator after the cleaning process. Whether in the isolator or on the BI, the organisms may not necessarily be in the form of a strict monolayer, but they must be presented such that the  $\text{VH}_2\text{O}_2$  can access the population with minimal restrictions. Therefore, it is important to select BIs that have been designed for  $\text{VH}_2\text{O}_2$  decontamination processes.

The authors of a recently published paper<sup>1</sup> demonstrated that BIs not optimized for a decontamination process can result in nearly every possible outcome when exposed to a  $\text{VH}_2\text{O}_2$  cycle. The authors obtained BIs from 4 manufacturers, with each

BI lot labeled for a purpose other than  $\text{VH}_2\text{O}_2$  decontamination (BIs varied in test organisms and spore counts). The BIs of each type were placed into the isolator and treated to a  $\text{VH}_2\text{O}_2$  process. The data presented in Table 1 is a summary of their results. (Note: The BIs in these experiments were not placed in the isolator such that they can be considered true replicates.) The data demonstrate results ranging from complete survivors with the BIs from manufacturer A to complete kill with BIs from manufacturer C and D, and fractional results in the remaining BI lot. These data demonstrate that it is crucial to use BIs that provide meaningful information about the process and that have been designed for use with  $\text{VH}_2\text{O}_2$ .

**Table 1. Summary of Data from Non-optimized BIs\* Exposed to a  $\text{VH}_2\text{O}_2$  Process<sup>1</sup>**

Manufacturer	A	B	C	D
BI's growth "+"/BI total	45/45	30/45	0/39	0/42


\*BIs that were designed for purposes other than  $\text{VH}_2\text{O}_2$

## $\text{VH}_2\text{O}_2$ BI Components

A BI is a system whose performance is dependent on the sum of its components. BIs properly designed for  $\text{VH}_2\text{O}_2$  consist of surrogate spore formers, the spore carrier, and the primary packaging if utilized. Each of these are discussed below.

### Spores

The surrogate organisms must be non-pathogenic and traceable to a recognized culture collection such as the American Type Culture Collection (ATCC). Spores of *Geobacillus stearothermophilus* (and to a lesser extent *Bacillus atrophaeus*) are recognized for use in  $\text{VH}_2\text{O}_2$  BIs as they have greater resistance to the process than the bioburden. There are two strains of *G. stearothermophilus* in use, ATCC #12980 and #7953. ATCC #12980 is more commonly used and provides a greater challenge than spores of the ATCC #7953 strain. Spores of *B. atrophaeus* (ATCC #9372) are utilized to a lesser degree and provide the lowest challenge to the process of the three strains discussed. Cleanliness of the spores is critical as cellular debris and growth media components can protect the spores. Figure 1 displays two states of cleanliness of a spore dilution. BIs made from dirty spore preparations may result in tailing or irregular performance.

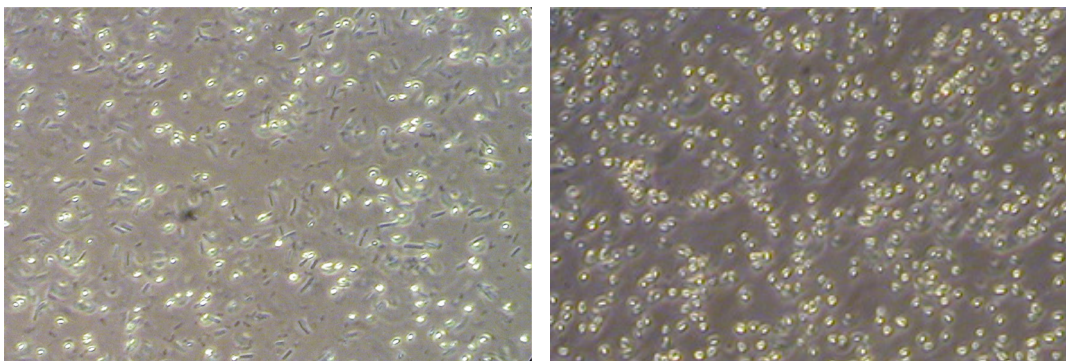

**Figure 1. Images of an Unclean Spore Preparation (left) and a Clean Spore Preparation (right), as viewed under a Phase Contrast Microscope**

## Spore Carrier

The spore carrier is simply the material on which the spores are deposited and is the vehicle in which the spores are delivered to their test locations in the isolator. A carrier material must not chemically react with or absorb the decontamination agent and must not impact spore viability. The majority of BIs produced globally for sterilization modalities utilize cellulose (i.e., paper filters) as the spore carrier, however for  $VH_2O_2$  processes, this material violates the above-mentioned criteria. The primary material in use today for  $VH_2O_2$  is stainless steel of various shapes (Figure 2.)

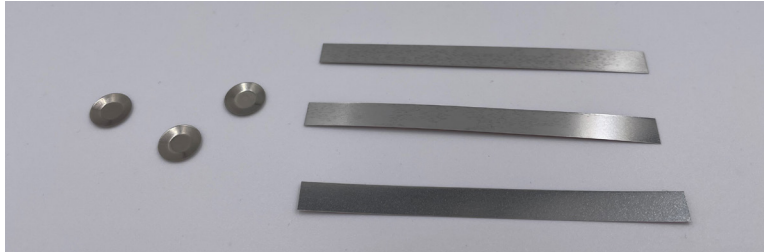


Figure 2. Stainless Steel Spore Carriers Typically Used in BIs for  $VH_2O_2$

The spore carrier should promote (to the greatest extent possible) the even distribution of spores across its surface. The ideal spore presentation is a complete monolayer; however, achieving this under the current recommendations ( $10^6$  spores placed over a small surface area) is not possible. The edge effect, which is the piling of spores on the outer boundary of the inoculated area (Figure 3.) will always be present and is generally not a concern, provided this feature is not exaggerated.

Although  $VH_2O_2$  may not penetrate well, it can certainly penetrate to some degree into smaller spore aggregates. If this were not true,  $VH_2O_2$  would never inactivate a BI and would not be suitable as a decontamination agent. The use of scanning electron microscopes (SEM) can provide great detail of spore distribution on a carrier, allowing the user to make some rudimentary and subjective predictions about the BI's performance, but it is not recommended to use this tool as part of a routine incoming inspection process. A study conducted at Mesa demonstrated the limitations of predicting BI performance based on SEM images<sup>2</sup>.

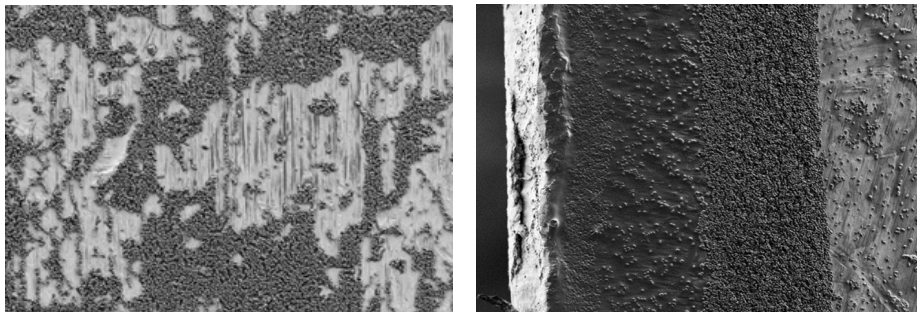


Figure 3. SEM Images of spores presented in a monolayer (left) and spore piling on the edge of the inoculum (right)

## Primary Packaging

The primary packaging contains the inoculated carrier and protects it from external contamination. The requirements for the packaging material are similar to those of the spore carrier in that it cannot chemically react with or absorb the decontamination agent and must not impact spore viability. Additionally, the packaging material must be permeable to  $\text{VH}_2\text{O}_2$ . The primary material in use today for  $\text{VH}_2\text{O}_2$  is Tyvek®, which is a commonly used packaging material for medical devices. It has been reported<sup>3</sup> that Tyvek does increase the resistance of the BI when compared to naked (unpackaged) versions of the same BI. These studies theorize that Tyvek allows the gas form of  $\text{VH}_2\text{O}_2$  to penetrate, while blocking the liquid form of the agent. For this reason, the use of naked BIs with fogging or misting systems may be more appropriate as this “vapor” tends to contain more liquid of various particle sizes.

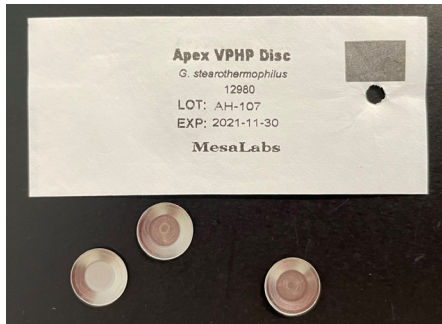


Figure 4. Image of spore discs and primary packaging

## Challenges with BIs in $\text{VH}_2\text{O}_2$ decontamination processes

To achieve meaningful information from the BIs and avoid unexpected BI results, certain steps should be considered including selection of the BI format and placement into the system. Important characteristics of BI components have been discussed above in detail; however, there are several options concerning BI format including packaged discs, unpackaged ribbons and Self-Contained BIs. A Spore News article<sup>4</sup> discusses the scenarios where one BI format may be advantageous over another, as well as techniques for proper placement in the isolators. If the BIs are not positioned properly, the spores may be shielded from the  $\text{VH}_2\text{O}_2$ , which increases the likelihood of a positive BI result.

Perhaps the main disadvantage of BIs as compared to other indicator systems is the time-to-results. To test for spore viability after a process, the spores on the BI must be cultured and incubated. The current regulatory expectation is an incubation period of 7 days, although a reduced incubation time is possible. In comparison, the results of a chemical indicator (CI) are known immediately post treatment, which makes them a valuable tool during the early stages of cycle development. However, despite the longer time-to-results, the BI is the only tool that can give a direct measurement of a spore log reduction (SLR) during the validation process as required by the regulatory agencies.

Manufacturers of BIs for  $\text{VH}_2\text{O}_2$  generally provide a “D-value” as a measure of resistance. By definition, a D-value is the time required to achieve one SLR under steady state conditions.  $\text{VH}_2\text{O}_2$  “D-values” are determined in a test isolator under stated conditions; however, these conditions are not at steady state for the duration of the process, nor do they account for all the variables that occur during the process. Agalloco<sup>5</sup> provides a detailed discussion on the physics of vapors (existing as both a gas and liquid) and the rationale for steady conditions not being met with  $\text{VH}_2\text{O}_2$  processes. Certainly, SLRs can be measured under the BI manufacturer’s stated conditions, but as the conditions of the test isolator change during the cycle, the killing rate (i.e., D-value) also changes. This is true in any  $\text{VH}_2\text{O}_2$  decontamination process. BI users are cautioned to not put too much weight on the manufacturer’s stated “D-value” for  $\text{VH}_2\text{O}_2$  decontamination processes, if for no other reason than the manufacturer’s test isolator will perform differently than the end user’s system. Therefore, the “D-value” presented by the manufacturer should only be considered a reference value.

### “Rogue” BIs

What exactly are “rogue” BIs, do they exist and if so, how prevalent are they in a BI lot? The term “rogue” BI was coined to describe an unexpected growth-positive BI unit that seemingly exhibits hyper-resistance to the process and is not representative of the BI lot. This growth positive BI unit is unexpected based on previously collected data in the same test locations often leading the practitioner to conclude that it is an outlier result. Unfortunately, it is not possible to directly prove whether a BI is “rogue” either before the treatment or afterwards when it results in spore outgrowth.

Unexpected positive BI results occasionally occur in traditional sterilization processes (moist heat, ethylene oxide, dry heat), but the term “rogue” BI hasn’t been used to describe these results. It seems the term was born out of the  $\text{VH}_2\text{O}_2$  decontamination process. “Spores don’t lie” is the term commonly used in traditional sterilization processes and simply means that if a BI is growth-positive after treatment (barring any laboratory error) then insufficient lethality was delivered during the process as demonstrated by spore outgrowth. But then, surface decontamination is not equivalent to terminal sterilization, so perhaps “rogue” BIs do exist for these processes.

The cause of “rogue” BIs has been attributed to very low levels of BIs within a lot that exhibit poor quality in the spore presentation on the carrier (i.e., extracellular debris, spore clumping or piling). The perfect BI lot would be one in which all spores on each carrier would exist in a perfect monolayer, which does not exist on BIs that contain a high spore density (e.g.,  $10^6$  spores on an 8mm disc). However, as noted above,  $\text{VH}_2\text{O}_2$  can penetrate smaller aggregate of spores as pictured in Figure 3, for if this were not true, then every BI in the lot would perform as a “rogue” BI.

Being that it is not possible to directly prove whether a BI is “rogue” either before or after the treatment, how does a user gain confidence on the suitability of the indicators prior to their use? BI lots can be evaluated as-a-whole either by the manufacturer (during the “D-value” assessment process) or by second- or third-party labs for general suitability in their process, including the presence of “rogue” BIs.

In a 2005 case study, Templeton and Hillebrand<sup>6</sup> describe in detail their approach in addressing the potential problems with unexpected positive BI results. In part, they performed a “rogue” detection test” on BIs from two suppliers. One lot showed a 1-2% “rogue” BI result while the other lot was less than 0.3%. What is an acceptable level? If the rate is 2% then perhaps the BI lot is unsuitable for use in this process, after all, can 2% of a BI lot really be considered “rogue” as they appear to represent a sizable portion of the batch. A result of <0.3% may be considered suitable for use with the knowledge that an occasional outlier is to be expected. The authors' approach of using multiple BIs at critical locations allowed them to demonstrate the necessary SLR even if/when an occasional unexpected positive BI resulted. Spore news Vol. 9 No. 4<sup>7</sup> provides additional details on the use of replicate BIs.

## Conclusions

Demonstrating the success of a  $VH_2O_2$  decontamination process is certainly more complex than traditional sterilization processes. This is due to poor agent penetration, non-steady state/multi-phasic conditions during treatment, and microorganism layering on isolator surfaces and BIs. Despite these challenges, proving a defined SLR value is not only possible, but it is also the current regulatory requirement. Additionally, the use of properly designed BIs will greatly increase the likelihood of a successful validation. Although there are no standards for the manufacture and performance of biological indicators for  $VH_2O_2$  decontamination processes, the details provided in this paper reflect proven and accepted practices. The biological indicator is the gold standard for measuring a spore-log-reduction as it is the only indicating system that actually contains viable spores.

## References

1. Macellaro, A; et al., Evaluation of Biological Indicator Spores as Tools for Assessment of Fumigation Decontamination Effectiveness. *Applied Biosafety*, 2015, 20 (4):183-191
2. Spore News Vol. 10, No. 6, SEM Analysis of VPHP Biological Indicators
3. Reich, R; Caputo, R. Vapor-Phase Hydrogen Peroxide Resistance of Environmental Isolates. *Pharmaceutical Technology*, Aug 2004: 50-58
4. Spore News Vol. 9, No. 5, End User---Proper BI Placement During VHP Decontamination Cycles
5. Agalloco, J; Vapor Phase Sterilization and Decontamination. *Handbook of Validation in Pharmaceutical Processes* 4th Edition, 2021 Chapter 21.
6. Templeton, P; Hillebrand J. Case Study: Isolator Sanitization Cycle Development, Validation and Revalidation. *Abstract of a paper presented at the ISPE Isolation Barrier Forums in Washington and Prague 2005*
7. Spore News Vol. 9, No. 4, Using Replicate BIs to Evaluate Biodecontamination Cycles in Isolators